

Rubber Fab Technologies Group's

CONTROLLED COMPRESSION GASKETS

Controlling the
Compressive Load to
Provide the Ultimate
Surface I.D.




Introducing the Controlled Compression Gasket

A Superior Hygienic Seal

The Controlled Compression Gasket cannot be over-tightened!

The CCG is composed of the elastomer of your choice with a ring of ultem® vulcanized inside the seal from the ring O.D. to the locating ring. This **non-compressible** polymer controls clamp compression to **exactly** the correct amount, allowing the properly positioned seal to make a perfect surface I.D. with the ferrules. **No intrusion of the hygienic seal** in the product flow is possible.

The Controlled Compression Gasket literally transforms the tri-clamp connection into a **perfect I.D. connection**. Because the ultem ring acts as a **positive mechanical compression stop** when clamping, the clamp will not loosen due to seal creep, aging or ΔT cycles. This is true for any **standard or high pressure clamp** regardless if the clamp is secured with a wing nut, standard nuts or Rubber Fab's Torque-Rite®

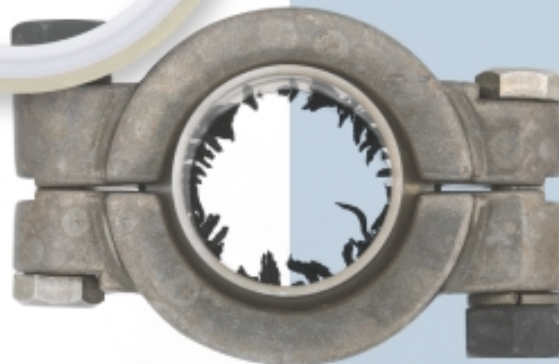

Over-tightening causes hygienic seal intrusion into the piping I.D. which can result in tearing and sticking resulting in system contamination.

Rubber Fab's Controlled Compression Gaskets act as a positive mechanical stop making "over-tightening" virtually impossible.

- Hygienic seal meets ASME-BPE intrusion requirement of $\pm .020$.
- Works with virtually any manufacturer's ferrules, fittings, clamps or combination thereof.
- No torque requirement. Tighten until it stops (we suggest wetting the gasket with isopropyl alcohol during installation to insure uniform compression, especially the 1" CCG).
- The CCG hygienic seal will allow for slight misalignment and ferrule face anomalies.
- Available in EPDM, FKM and silicone.
- Can be tabbed for traceability or non-tabbed with laser engraved batch and lot information.
- FDA compliant, 3A and Class VI tested.
- First major gasket design patent issued in the last 60 years.



Controlled Compression Gaskets

- Hygienic seal material will not extrude into I.D. maintaining perfect surface I.D.
- Transforms tri-clamp fittings into a perfect I.D. connection.
- Maintains seal after thermal cycling.
- No re-tightening required.
- Complete control over amount of compression load on seal material.
- Creates a static and leak tight seal.
- Makes over-torquing impossible.
- Superior cleaning, draining and sanitization guaranteed.
- Certain elastomers can withstand up to 100 steam cycles.
- Prevents soil retention and fluid hold-up.
- Tabs indicate lot, batch, manufacture date, size and materials used.
- Constant lot and batch traceability.
- Tabs insure no potential human contamination during installation.
- Works with standard and high pressure clamps.
- 1" to 3" sizes available (excluding 2-1/2").
- Complete traceability with color coded system.
- Adjusts for misalignment.
- Available standard or tabbed.
- Metal Detectable Elastomers available.

© Torque-Rite and ADI Free are registered trademarks of Rubber Fab Technologies Group.

Controlled Compression Gasket: United States Patent Number 6,857,638.

Torque-Rite: United States Patent Number 6,082,941.

Rubber Fab Technologies Group's Metal Detectable products are patent pending.

© Ultem is a registered trademark of GE.



± † ¶ ‡ ■ ○
Rubber Fab Technologies Group
Limecrest Road

P.O. Box 626 • Andover, NJ 07821
phone: 973-579-2959 fax: 973-579-7275

www.rubberfab.com

Perfect...Every Time


Because there is **no unsupported hygienic seal intruding into the product flow**, the seal's **service life is greatly extended**. With the minimum surface contact area exposed, particulates, TOCs and other extractables are dramatically reduced.

- Compression of the hygienic seal surface is more than adequate to compensate for imperfections in ferrule faces. The ultem® ring is ductile enough to allow for piping misalignment.
- The lettering on the ultem ring indicates the elastomer type. Ring can be a sanitary Type I standard style or tabbed. Tabbed ring provides permanent record of batch and lot number as well as date of manufacture for tracking. The tabbed ring allows customer to eliminate human contact with the hygienic seal surface and an area to record date of installation on the reverse side of the tab.
- Ultem is an FDA compliant, USP Class VI approved, non

compressible polymer. Maximum operating temperature rating is 450°F. Controlled Compression Gaskets can be autoclaved and depending on the elastomer selection can withstand up to 100 steam cycles.

Meeting the Stringent Standards For Purity

- U.S. Pharmacopeia Class VI Certification
- Cytotoxicity Criteria
- Title 21 CFR 177.2600
- USDA and 3-A Sanitary Standards
- Current Good Manufacturing Practices (CGMP)
- ASME-BPE Standards
- Animal Derived Ingredient Free

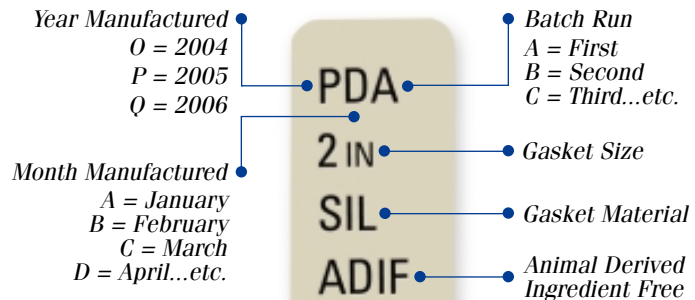

Creating a Sanitary Tracking and Coding System

Rubber Fab Technologies Group is a member of:

Distributed by:

